

CAGEWASH 3

ALKALINE DETERGENT CLEANER

Concentrated Alkaline Detergent Cleaner

Heavy duty alkali detergent ideal for use on cage washers, bottle washers and rack washers for the removal of organic and greasy soiling from caging and associated equipment.

Economical and versatile with multi-surface use, suitable on both hard and soft water applications. Has excellent rinsing capabilities and effectively removes heavy, baked on organic and greasy deposits often seen on larger caging, and racks.

Suitable for use on low temperature washing below 50°C.

Application:

1. Place feed tube into bottom of container and check automatic feed equipment is switched on.
2. Inject at 4ml per 1 litre of water at a working temperature of 45 – 60 °C.

Material Compatibility:

Cagewash 3 is compatible with the following materials:

Stainless steel, aluminium, anodised aluminium, chromium plated material, synthetic materials, rubber, latex, glass, ceramics, copper, brass, and metal alloys.

Warnings & Cautions:

- Do not add any other chemicals, such as bleach or detergents as this may destroy the effectiveness of this product
- Wear goggles, plastic apron and rubber gloves when changing containers.
- Consult the manufacturers recommended reprocessing instructions to determine cleaning agent compatibility and any additional procedures necessary to clean their particular medical device
- Store in a cool place between 5 and 25°C
- Dispose in accordance with local authority guidelines



Available in 5L, 10L, 20L, and 200L